

**RODENT EUTHANASIA BY CO₂ INHALATION
IS PERFORMED BY USING GRADUALLY
INCREASING CONCENTRATIONS OF CO₂
AT RATES INDICATED BELOW:**

Species	Cage Type	Flow Rate L/min
Mice	Static (small)	2
	Static (large)	6
	IVC (w/o lid)	3
	IVC (w/ lid)	4
Rat	Static	8
	IVC (w/o lid)	11
	IVC (w/ lid)	13

SC #089.3
Eff. 3/17

4x6

**Finish size- add 1/2" border around all 4 sides of table
Print and laminate to be displayed at all CO₂ stations**